

COLIFORM WATER SAMPLING & ANALYSIS VIALS

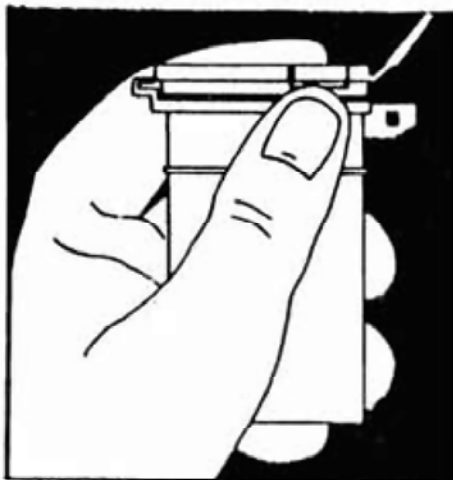


FIG. 1

To open, firmly hold container and apply enough pressure against thumb tab to break pre-use, arrow locking sterility seal.



FIG. 2

Grab Lock-N-Seal Latch and move lid away from collection stream. Exercise care not to touch inside of lid or vial.



FIG. 3

Collect sample using proper sample collection techniques.

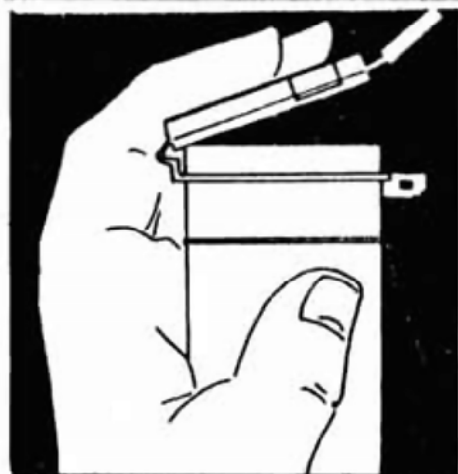


FIG. 4

To close, press down on lid from behind using forefingers. Snap shut and firmly press to sealing ring.



FIG. 5

Snap down Lock-N-Seal Latch to avoid accidental opening during transport.



FIG. 6

Pull custody tie thru hole in latch & complete information on sample label. (Custody tie & pre-applied labels are optional.)